






MDF-DU901VHA-PA

VIP°ECO

Large Volume -86°C Ultra-Low Temperature Freezer with Natural Refrigerants

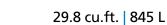

MDF-DU901VHA-PA, 115V









The MDF-DU901VHA-PA is a large volume high performance ultra-low temperature freezer ideal for use in biorepositories or facilities where bulk storage of sensitive biologicals are managed. This freezer delivers optimum performance over a range of setpoints from -50°C to -86°C, adjustable in 1°C increments. It operates with natural refrigerants at 115V power and can pass through a standard 36" wide doorway for new or replacement installation; (may require temporary removal of exterior stand-off brackets).

Natural Refrigerants, Intelligent Compressors

The robust refrigeration system is powered by a new generation of high performance inverter compressors with natural refrigerants. These fast temperature recovery following ambient conditions.

compressors are supervised by proprietary algorithms in the main controller firmware to maintain the most efficient energy consumption. Reserve refrigeration power assures door openings and tolerance for high

Microprocessor Controller

An integral microprocessor controller combines setpoint security, monitoring, predictive performance and data logging functions into a single eye-level display. The large color LCD touchscreen can be controlled with a gloved hand. A USB port permits downloading of operating history as desired.

Energy Efficient Performance

The MDF-DU901VHA-PA combines upright freezer cabinet design, vacuum composite insulation and variable speed compressors that operate independently and only on demand. All natural refrigerants support facility sustainability objectives without compromising reliability, performance or reserve power needed to tolerate real world conditions. Operating costs are minimized while stored product protection is maintained.



Large Volume Efficiency

As sample volumes in biorepositories and biobanks increase exponentially, the high capacity storage chamber offers space for expansion and overflow from crowded smaller freezers.

High Capacity Storage, **Minimal Footprint**

The upright cabinet design includes a patented VIP Plus insulation composite which creates maximum interior space efficiency. It is designed and tested to tolerate high ambient temperatures in warm labs or hallways. The upright cabinet orientation holds up to 672, 2" boxes or 448, 3" boxes within a 10.79 sq.ft. (~1.0 sq.m) footprint.



Ergonomic Interface

The EZlatch door handle (patent pending) is central to the inventory access system. It permits secure, onehand opening and closing with a minimum of effort while preventing frost build-up. Positive latching insulated inner doors supplement outer door insulation. A manual vacuum relief port permits quick re-entry.




VIP°ECO -86°C Ultra-Low Temperature Freezer with Natural Refrigerants



VIP Plus Insulated Panels

Inside-out engineering starts with a VIP Plus insulation composite which results in an efficient, thin-wall cabinet. Insulated inner doors improve temperature uniformity and cold air loss during door opening and extend warm-up times during power outages.


Temperature Uniformity Data:

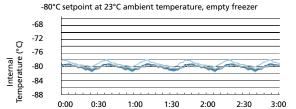

Time (hours:minutes)

Dimensions

Unit: inch | mm

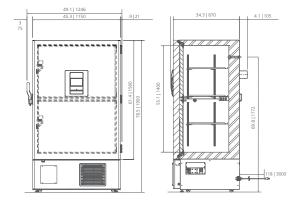



PHC Corporation of North America 1300 Michael Drive, Suite A, Wood Dale, IL 60191 Toll Free USA (800) 858-8442, Fax (630) 238-0074 www.phchd.com/us/biomedical

er		
Model Number		MDF-DU901VHA-PA
External Dimensions (W \times D \times H) 1)	inches mm	45.3 × 34.3 × 78.5 1150 × 870 × 1993
Internal Dimensions (W \times D \times H)	inches mm	39.8 × 23.6 × 55.1 1010 × 600 × 1400
Volume	cu.ft. liters	29.8 845
Net Weight	lbs kg	723 328
Capacity (2 * boxes)	qty	672
Capacity (3 * boxes)	qty	448
Inventory Racks (2" boxes)	qty	28
Inventory Racks (3" boxes)	qty	28
Performance		
Warranty 2)		5 years parts and labor
Cooling Performance 3)	°C	-86
Temperature Setting Range 3)	°C	-50 to -90 in 1° increments
Temperature Control Range 3)	°C	-50 to -86 in 1° increments
Control	C	30 to do iii i iideinens
Controller		Microprocessor, touchscreen data entry, password protected
Display		LCD color touchscreen
Temperature Sensor		Pt-1000
Refrigeration		
Refrigeration System		Synchronized variable speed cascade system
Compressors	W	(2) 1000, variable speed
Refrigerant		High-stage R-290 natural refrigerant
		Low-stage R-170 natural refrigerant 3.1 78, rigid polyurethane foam (PUF)
Insulation Thickness, Material	inches I mm	+ VIP Plus vacuum insulated panels
Construction		
Exterior Material		Painted electrogalvanized steel
Interior Material		Powder coated electrogalvanized steel
Outer Door	qty	1 (key lock with provision for optional padlock)
Inner Doors	qty	2 (insulated, ABS, stainless steel frames, positive latches)
Shelves	qty	3 (adjustable)
Shelf Dimensions (W × D)	inches mm	38.9 × 20.9 988 × 533
Max. Load per Shelf	lbs kg	110 50
Max. Load - Total Freezer		1266 414.4
	lbs kg	1: lower left side
Manual Vacuum Relief Access Ports	qty	
	qty	2: back wall and chamber floor
Access Port Diameter	inches mm	0.6 17
Casters	qty	4 (2 leveling feet on front base)
Alarms	(V =	Visual Alarm, B = Buzzer Alarm, R = Remote Alarm)
Power Failure		V-B-R
Temperature (High and Low)		V-B-R
Sensor Failure		V-B-R
Door Open		V-B
Battery Check		V
Remote Alarm Contacts		DC 30V 2A (normally open, normally closed, common)
Electrical and Noise Level		
Power Supply	(-) (h	115V, 1Ø, NEMA 5-20P, requires NEMA 5-20R receptacle
Noise Level 4)	dB(A)	52
Options		
<u> </u>		AIDE CIT SILL
Inner Door Kit with 4 Smaller Doors 5)		MDF-9ID-PW
Liquid CO₂ Back-Up System		MDF-UB7-PW
Liquid Nitrogen Back-Up System		MDF-UB7-PW-LN2
Temperature Recorder		
Circular Type	6", 7 day circular	MTRC954
Chart Paper	52 charts per box	C7100386REV
Ink Pen	pack of 6	R252
Optional Communication System		
Wireless, Cloud-Based,		LIAL (SAL TO A CO.
Automatic Data Management		LabAlert® Monitoring System

- ¹⁾ Exterior dimensions of cabinet excluding handle, rear stand-off brackets and other external projections. Consult product web page for doorway entry instructions, less than 36": www.phchd.com/us/biomedical/preservation/ultra-low-freezers/mdf-du901vhapa
- 2) Current warranty offered at time of printing and may be subject to change; US and Canada only 3) Air temperature measured at freezer center, ambient temperature +30°C, no load
- 4) Nominal value, background noise 20 dB(A)
- 5) Installation of small inner door kit may affect usable storage capacity

Specifications are subject to change without notice.

For latest specification information contact PHC Corporation of North America at info@us.phchd.com. Performance data herein is based on independent testing at time of publication.

